λ /HindIII DNA for long ssDNA preparation kit

Cat. #	Product	Size
DS670	λ /HindIII DNA for long ssDNA preparation kit	600 µl


Storage Conditions: This product should be stored at -80°C.

Description:

This product is the component of Long ssDNA Preparation Kit for 10kb (Code #DS635). The approximate size of ssDNA can be estimated by this product.

Product Usage:

- 1. Aliquot 20 µl to new tube.
- 2. Heat at 70°C for 5 min.
- 3. Chill on ice for 1 min.^{*1}
- 4. Subject to agarose gel electrophoresis with 1×TAE. *2, *3
 - *1 After heating and chilling, quickly load into wells of the gel and promptly start electrophoresis.
 - *2 The mobility of ssDNA is more changeable depending on the sequence than dsDNA, therefore the bands indicates rough size.
 - *3 23130 bases and 6557 bases may appear as double-band.

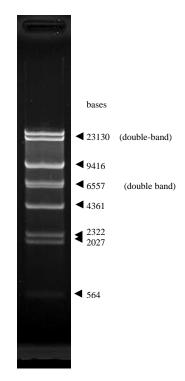

1.2 % agarose gel in $1 \times TAE$

1

Related Products:

DS635	Long ssDNA Preparation Kit for 10kb
DS650	Long ssDNA Gel Extraction Kit for 10kb

ver. 1.0